

RESULTING A RAPID STREP SCREEN OR RSV AG IN CERNER



1. Open Microbiology Result Entry from the Appbar.
2. Enter the accession number into the Entry box or scan the barcode.

Accession: Proc

3. The Microbiology Result entry screen will open.

4. In the Entry box, enter LOTQC

Entry:

Biochemical	Coded Result	Coded Result Desc	Freetext Result	Status	Date/Time	User
Lot#				Pending		E233341
Exp Date				Pending		E233341
Pos Control				Pending		E233341
Neg Control				Pending		E233341

OK Cancel Mark Unmark

The Add Biochemical –LotQC box will pop up.

Free text or scan a barcode to result the Lot# and Expiration Date.

Biochemical	Coded Result	Coded Result Desc	Freetext Result	St
Lot#			123456	Pendir
Exp Date			01/01/2020	Pendir
Pos Control	A	Acceptable		Pendir
Neg Control	A	Acceptable		Pendir

Perform Verify

and then use the pull down arrows to choose “A” for Acceptable for both the Pos and Neg Controls. This represents the original QC that was done on this kit.

The result for QC will now show up on the bottom half of the screen.

This is not a visible result to the doctors.

Update Dt/Tm	Tack ID	Status	#	Enter	Lot(s)	Biochemical Group	Res	Test	Lot Number
12/17/2015 14:24	E233341	Verified		Lot#		Test: 123456	Res:	Test: 01/01/2020	Expiration Date
12/17/2015 14:24	E233341	Verified		Exp Date		Res: A	Test: Acceptable	Positive Contro	
12/17/2015 14:24	E233341	Verified		Pos Control		Res: A	Test: Acceptable	Negative Contro	

5. Now go back to the Entry box and type in the word “FINAL” and the Final Report screen will pop up.

Final Report

Page:

Get Up Get Down Get Left Get Right Print Out


Print Out Deep Edit Erase Select Row

Chat View

Route Cancel Route Perform Verify


6. The following responses must be used to result in the result box:

TEST	ORDER	RESULTS	MEANING
STREP SCREEN	GROUP A STREP AG	SAAN2	NEG -PRESUMPTIVE...
		SAAP	POS
		IPCA	INTERNAL QC OK
RSV	RSV ANTIGEN DETECTION	NRSVIA	NEG
		RSVP	POS
		IPCA	INTERNAL QC OK

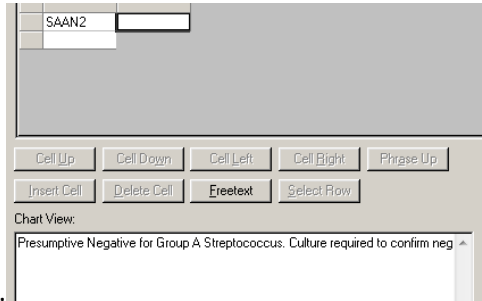


Choose the appropriate result, for example...

And press enter. Press enter again to go to the box below. This is where the code IPCA must

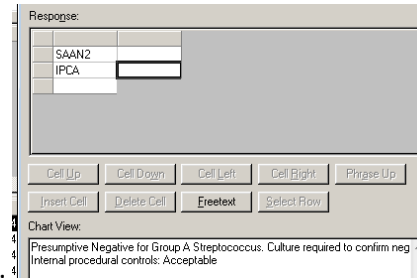


always be entered.

At this point, check the box for the Chart View and make sure it says exactly what you want to report. This is what will be visible to the doctors. If this is correct,

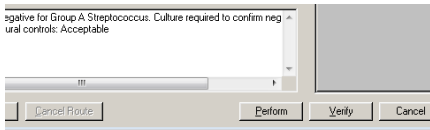

choose Verify.

7. If the test result was negative for the rapid strep screen, then a Throat Culture will automatically be ordered on the same accession number. Note: This will not work properly if Perform is used and then later Verify is pressed. All that needs to be done now is to reprint the barcode to send it out.

8. To Reprint a label, choose the Reprint Label app , scan or enter the accession number, and a new label will print showing both the Throat culture and the rapid strep screen.